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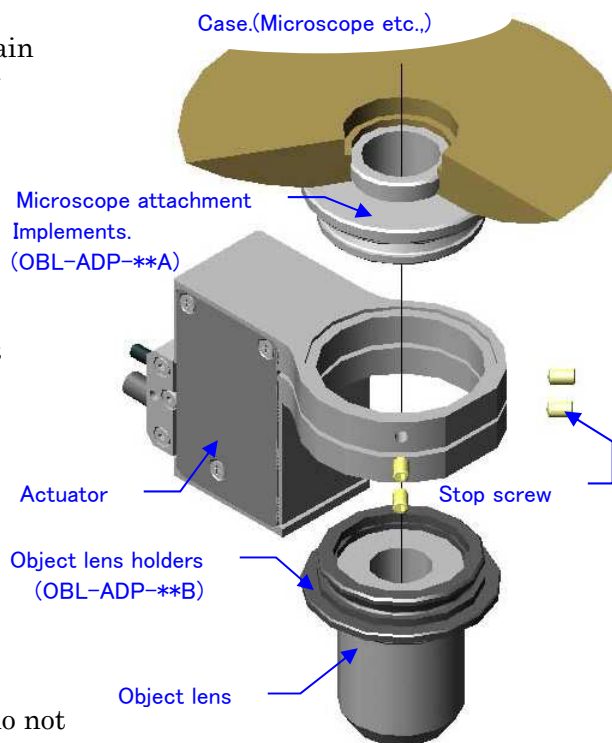
SIGMA FINE STAGE (The Actuator for Object Lenses)

- SFS-OBL-1/-2 - Installation manual


I. Installation

- (1) Take out the actuator from the storage case and refrain for use the storage case for storing and transporting the actuator.
- (2) An object lens is certainly thrust into an object lens holder, and it fixes.
- (3) A microscope attachment implement is certainly thrust into a case (microscope etc.), and it fixes.
- (4) An object lens holder is inserted in a main part, two stop screws are tightened, and it certainly fixes.
- (5) The actuator is inserted in a microscope attachment implement, two stop screws are tightened, and it certainly fixes.

※ Precautions for Use

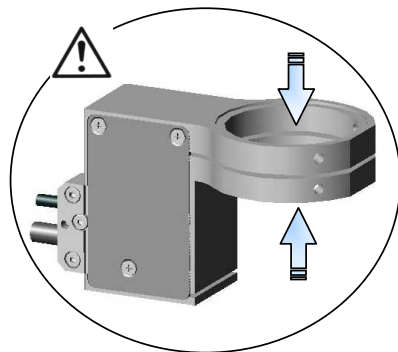- In case you treat the actuator, be careful enough not to give an impact to the actuator or hit it. Otherwise, the displacement mechanism may be damaged. Do not give pressure to the fixed part and movable part of the actuator.
- In case you operate a case (microscope etc.), please do not give an impact to the tip of an object lens.

⚠ Do not touch during operations.

⚠ Should be rounded a case(microscope..) properly in your environment.

⚠ Always check that the power to the controller is turned off before connecting. If the power is turned on, the actuator may not operate correctly or become damaged.

II. Environment of Use

Use the stage in the following environment.

Temperature 10°C~30°C

Humidity 20%~60% (With no dew condensation)

REMARKS: The metal tends to elongate as temperature rise. For more accurate positioning, minimize the temperature change as much as possible during usage.

(Recommended temperature $\pm 1^\circ\text{C}$)

III. Storage

When storing, fix the actuator to the special case using screws. Do not give an impact to the stage at this time. Store the stage in the following environment.

Temperature : 0°C~40°C

Humidity : 10%~85% (With no dew condensation)